

M-011 Linear Slide with DC/Piezo Drive Compact, Nanometer-Precise Piezo Drive, High-Accuracy Guiding



M-011.D01 translation stage

- Travel Range to 15 mm
- Compact Side Drive
- Straightness/Flatness ≤0.2 µm
- PZT Drive for Scanning and Tracking Applications
- 0.1 µm Resolution with Closed-Loop DC Motor
- 5 nm Resolution with Closed-Loop PZT Drive

© Physik Instrumente (PI) GmbH & Co. KG 2008. Subject to change without notice. All data are superseded by any new release. The newest release for data sheets is available for download at www.pi.ws. R1 10/03/10.0

M-011 ultra-high-precision magnetically-coupled stages use the force of integrated magnets to preload the bearing. This magnetic preload results in extremely uniform and smooth motion with minimum friction. Unlike conventional stages, where two bearings with limited parallelism guide the carriage (inducing runout and rotational errors) in M-011 stages, only one of the two linear bearings has a guiding function (V-groove) while the second bearing is for support only (U-groove).

Ultra-Straight Motion

The coupling between the stage and the space-saving side-drive units (DC-motor drives, PZTs, micrometer drives) is not rigid but via mag-

Application Examples

- Microscopy
- Quality control
- Metrology

nets. This design allows only on-axis forces (drive direction) to be transmitted to the stage; torque-induced positioning errors resulting from non-parallelism between the drive axis and the guiding axis are eliminated.

Six Different Versions

The basic version, the M-011.00, is equipped with a precision manual micrometer providing a sensitivity of 1 μ m. M-011.D01 is equipped with a closed-loop DC-motor drive providing 0.1 μ m minimum incremental motion.

High-Resolution Piezo Option

For both the manual and motorized version, closed-loop and open-loop piezo drives are available. They provide 5 nm minimum incremental motion over a travel range of 30 μ m and allow for dynamic operation such as scanning and tracking. The closed-loop piezo drive provides repeatability of 60 nm (see the "Piezo Actua-

tors" section for further details on piezo actuators and recommended controllers). All stages can be cross stacked and combined with the M-052 Z-axis mounting bracket to provide multi-axis motion (vertical use of the M-011 only permitted with loads less than 0.1 kg, no extended Z-axis bracket for motorized Z-stage available).

Upgrades

M-011 stages without piezo drives or DC-motor drives can be upgraded at a later date (see ordering information).

Notes

See "Accessories" (p. 4-90 *ff*) for adapters, brackets, etc.

Ordering Information

M-011.00 Translation Stage, 15 mm

M-011.P0 Translation Stage, 15 mm, Manual

+ PZT Drive

Translation Stage, 15 mm, Manual + Closed-Loop PZT Drive

M-011.D01 Translation Stage, 10 mm, DC-Motor Drive

M-011.DP1 Translation Stage, 10 mm, DC-Motor + PZT Drive

M-011.DS1 Translation Stage, 10 mm, DC-Motor + Closed-Loop PZT Drive

Upgrades

M-011.U0 Upgrade Kit with Open-Loop PZT Drive

M-011.US Upgrade Kit with Closed-Loop PZT Drive

M-011.UD1 Upgrade Kit with DC-Motor Drive and Limit Switches (Factory Installed)

Ask about custom designs!



XY combination of two M-011.00 stages



M-011.00 translation stage









Technical Data

Model	M-011.00	M-011.P0	M-011.PS	M-011.D01	M-011.DP1	M-011.DS1	Units
Travel range	15	15	15	10	10	10	mm
Piezo fine travel range	-	30	30	-	30	30	μm
Min. incremental motion (piezo drive)	-	0.005	0.005	-	0.005	0.005	μm
Repeatability (piezo drive)	-	-	0.06	-	-	0.06	μm
Design resolution (DC Motor)	-	-	_	0.0035	0.0035	0.0035	μm
Min. incremental motion	1	1	1	0.1	0.1	0.1	μm
Unidirectional repeatability	-	-	-	0.1	0.1	0.1	μm
Bidirectional repeatability	-	-	-	2	2	2	μm
Backlash	-	-	-	2*	2*	2*	μm
Straightness / flatness per 5 mm	0.1	0.1	0.1	0.1	0.1	0.1	μm
Straightness / flatness full travel	0.2	0.2	0.2	0.2	0.2	0.2	μm
Max. velocity (motor)	-	-	-	0.7	0.7	0.7	mm/sec
Max. normal load capacity	1	1	1	1	1	1	kg
Max. push/pull force	7 / 5	7 / 5	7 / 5	7 / 5	7 / 5	7 / 5	Ν
Max. lateral force	5	5	5	5	5	5	N
Drive	M-623	M-623	M-623	M-227.10	M-227.10	M-227.10	
Piezo drive	-	P-840.20	P-841.20	-	P-840.20	P-841.20	
Encoder resolution	-	-	-	2048	2048	2048	counts/rev.
Drivescrew pitch	0.5	0.5	0.5	0.5	0.5	0.5	mm/rev.
Gear ratio	-	-	-	69.12:1	69.12:1	69.12:1	
Nominal motor power	-	-	-	2	2	2	W
Motor voltage	-	-	-	12	12	12	V
Mass	0.35	0.4	0.42	0.55	0.6	0.62	kg
Body material	AI, St	Al, St	Al, St	AI, St	Al, St	Al, St	
Recommended piezo controller	-	E-660, E-610	E-610	-	E-660 (p. 2-119), E-610	E-610 (p. 2-110)	
		E-500 System	E-500 System		E-500 System	E-500 System	(p. 2-142)
Recommended motor controller	-	-	-	C-843, C-848, C-863	C-843, C-848, C-863	C-843, C-848 (C-863 (p. 4-11)	p. 4-120 <i>ff</i>), 4)

Linear Actuators & Motors

Nanopositioning/Piezoelectrics

Nanometrology

Micropositioning

Hexapod 6-Axis Systems / Parallel Kinematics

Linear Stages

Translation (X) Vertical (Y) Multi-Axis

Rotary & Tilt Stages

Accessories

Servo & Stepper Motor Controllers	
Single-Channel	
Hybrid	
Multi-Channel	

Micropositioning Fundamentals

```
Index
```

*gearhead